





# Instructions for loading buckets

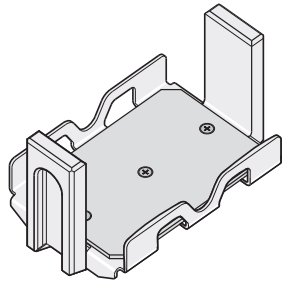


**5804 743.008 / 022638556**  
Set of 2  
DWP buckets for swing-bucket rotor A-2-DWP  
**5804 740.009 / 022638564**

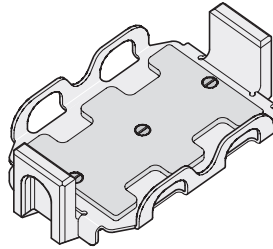


**5427 722.009 / 022634420**  
Set of 2  
MTP buckets for swing-bucket rotor A-2-MTP  
**5427 700.005 / 022634403**

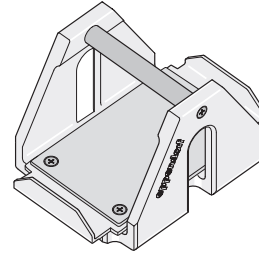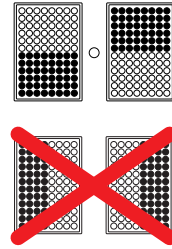

**5810 702.003 / 022638068**  
Set of 4  
MTP buckets for swing-bucket rotor A-4-62-MTP  
**5810 711.002 / 022638041**

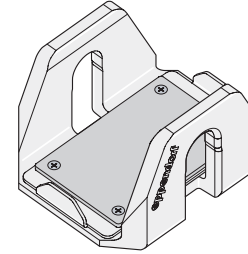


**5810 741.009 / 022638840**  
Set of 4  
MTP/Flex buckets for swing-bucket rotor A-4-81-MTP/Flex  
**5810 725.003 / 022638807**

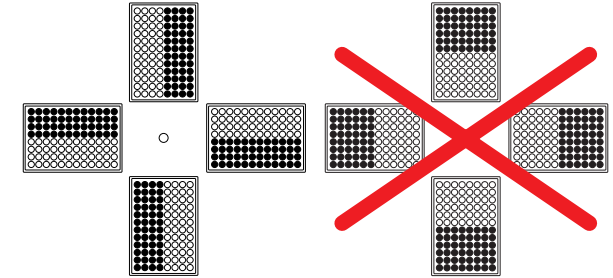


## Incorrect loading


- Unsuitable format, e.g. racks **too small**,

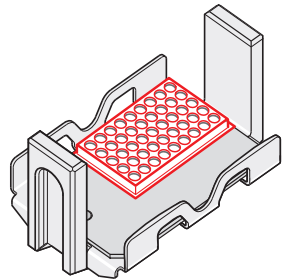

- microtiter plates **too wide**.

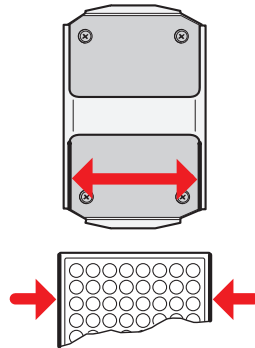


- Stacked microtiter plates **from several different** manufacturers,

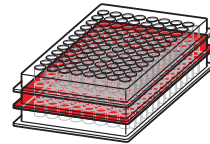

- **mixed loading** of deepwell plate and microtiter plates.



- Microtiter plates or deepwell plates **not inserted correctly**.

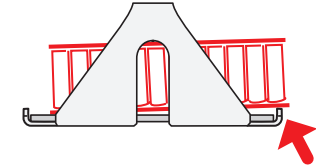

Deepwell plate buckets of swing-bucket rotor A-2-DWP, A-4-62-MTP and A-4-81-MTP/Flex may be loaded with plates (commercially available microtiter plates, culture plates and deepwell plates only), which do not exceed a weight of 380 g. The maximum load for the buckets of the swing bucket rotor A-2-MTP is 170 g. When using plates other than those named above, the user must first consult the Eppendorf Application Hotline. A manual loading and swing test are to be carried out

prior to centrifugation of plates or plate stacks that have not yet been used.

Maximum permissible load height:	
DWP bucket (A-2-DWP)	89 mm
MTP bucket (A-2-MTP)	29 mm
MTP bucket (A-4-62-MTP)	53 mm
MTP/Flex bucket (A-4-81-MTP/Flex)	60 mm

**eppendorf**  
*In touch with life*

Your local distributor: [www.eppendorf.com/worldwide](http://www.eppendorf.com/worldwide)

Eppendorf AG · Germany · Tel. +49 40 538 01-0 · E-Mail: [eppendorf@eppendorf.com](mailto:eppendorf@eppendorf.com) · Eppendorf North America, Inc. · USA · Toll free phone 800 645 3050 · E-Mail: [info@eppendorf.com](mailto:info@eppendorf.com)

**Application Support** Europe, International: [support@eppendorf.com](mailto:support@eppendorf.com) · North America: [support\\_NA@eppendorf.com](mailto:support_NA@eppendorf.com) · Asia, Pacific: [support\\_Asia@eppendorf.com](mailto:support_Asia@eppendorf.com)